

MACHEREY-NAGEL

QUANTOFIX[®] Relax

Automated reader for test strips

Rapid Tests



Perfect vision · Exact results

- Highest precision
- Objective readings
- Easy result administration



MACHEREY-NAGEL

www.mn-net.com



QUANTOFIX® Relax

A revolution in test strip analysis

The QUANTOFIX® Relax is a revolutionary step in test strip analysis. With this new system, MACHEREY-NAGEL offers an entirely objective method for reading QUANTOFIX® test strips, making the system particularly suitable if results are critical to your decision making. The system simplifies testing procedures and the decision processes by eliminating the subjectivity of visual readings and by providing higher accuracy. Especially if you test large numbers of strips, the QUANTOFIX® Relax helps and supports you with easy data administration and quick testing procedures. Additionally, the QUANTOFIX® Relax allows you to get actual quantitative results – no gradation and estimation of in-between values is necessary.

The QUANTOFIX® Relax can change the way you work and test. The system enables you to focus on your core processes and is the perfect tool for your testing procedures.

Speed up your analysis and save time with up to 50 analyses per hour!

Highest precision

- Quantitative results of complete measuring range
- Higher accuracy and no in-between value guesswork
- Assure your decision making

Objective readings

- No influence of external light or human color perception
- Reproducible results independent of the operator
- Increase safety and comfort

Easy result administration

- Immediate printouts
- Data can be transferred to a computer
- Save time to focus on your core business




Optimized usability



Auto start

Select the test, dip the strip and place it on the instrument. Without pressing any further button, the instrument starts the measurement automatically. Thus, operating is clean and easy.

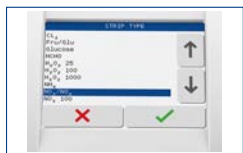

Graphic display

The graphic display shows all relevant information at a glance. The result is given with a test specific confidential interval. Thus, you can easily decide which further steps are needed for your process.



Touch screen

The touch screen makes using the instrument a real pleasure. All functions are easily accessed with a tap on the screen. You can also enter sample IDs with the on-screen keypad or an external keyboard. This makes documentation of results safe and easy.



Interface options

The instrument features an USB as well as a RS 232 interface. It can easily be connected to existing laboratory information systems and also to PCs.




Printer

The result is printed using a fast and silent internal printer. Thus, you can easily staple the results to production files for later check or QC procedures.



Easily change tests

With just two taps on the screen you can quickly change between your favorite tests. Select five different tests as your favorites out of a large variety of tests programmed into the instrument. All information is already programmed so you can immediately start the measurement when needed.

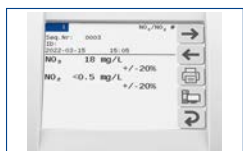

Memory for 200 measurements

Up to 200 measurements can be stored. Using internal functions, positive readings are easily filtered out, enabling further diagnosis whenever necessary.



Test strips for QUANTOFIX® Relax

The following tests are calibrated on the instrument and can be run with the instrument's software.

Test	Range	REF
■ QUANTOFIX® Ammonium*	10 – 350 mg/L NH_4^+	91315
■ QUANTOFIX® Ascorbic acid	25 – 1000 mg/L vitamin C	91314
■ QUANTOFIX® Chlorine Sensitive	0.1 – 10 mg/L Cl_2	91339
■ QUANTOFIX® Chromate	3 – 100 mg/L CrO_4^{2-}	91301
■ QUANTOFIX® Formaldehyde*	10 – 200 mg/L HCHO	91328
■ QUANTOFIX® Total iron	5 – 500 mg/L $\text{Fe}^{2+} \text{ } ^{3+}$	91330
■ QUANTOFIX® Total acid	2 – 5 g/L citric acid	91353
■ QUANTOFIX® Total sugar	55 – 700 mg/L fructose / glucose	91352
■ QUANTOFIX® Glucose	50 – 2000 mg/L glucose	91348
■ QUANTOFIX® Copper	10 – 300 mg/L $\text{Cu}^{+} \text{ } ^{2+}$	91304
■ QUANTOFIX® Nickel	5 – 100 mg/L Ni^{2+}	91305
■ QUANTOFIX® Nitrate 100	3 – 100 mg/L NO_3^- / 0.5 – 50 mg/L NO_2^-	91351
■ QUANTOFIX® Nitrate / Nitrite	10 – 500 mg/L NO_3^- / 0.5 – 80 mg/L NO_2^-	91313
■ QUANTOFIX® Nitrite	0.5 – 80 mg/L NO_2^-	91311
■ QUANTOFIX® Nitrite 3000	0.1 – 3.0 g/L NO_2^-	91322
■ QUANTOFIX® Peracetic acid 50	5 – 50 mg/L peracetic acid	91340
■ QUANTOFIX® Peracetic acid 500	50 – 500 mg/L peracetic acid	91341
■ QUANTOFIX® Peracetic acid 2000	500 – 2000 mg/L peracetic acid	91342
■ QUANTOFIX® Peroxide 25	0.5 – 25 mg/L H_2O_2	91319
■ QUANTOFIX® Peroxide 100	1 – 100 mg/L H_2O_2	91312
■ QUANTOFIX® Peroxide 1000	50 – 1000 mg/L H_2O_2	91333
■ QUANTOFIX® Phosphate*	3 – 80 mg/L PO_4^{3-}	91320
■ QUANTOFIX® Sulfite	10 – 500 mg/L SO_3^{2-}	91306
■ QUANTOFIX® Zinc	5 – 40 mg/L Zn^{2+}	91310
■ pH-Fix 0 – 14**	pH 1.0 – 13.0	92110
■ pH-Fix 0.0 – 6.0	pH 0.5 – 6.0	92115
■ pH-Fix 2.0 – 9.0	pH 2.0 – 9.0	92118
■ pH-Fix 3.6 – 6.1**	pH 3.6 – 6.1	92130
■ pH-Fix 4.5 – 10.0	pH 4.5 – 10.0	92120
■ pH-Fix 6.0 – 10.0	pH 6.0 – 10.0	92122
■ pH-Fix 6.0 – 7.7	pH 6.0 – 7.7	92150
■ pH-Fix 7.0 – 14.0	pH 7.0 – 13.5	92125

Together with the QUANTOFIX® Relax, test strips may not be used for any medical application.

* This product contains harmful substances which must be specially labeled as hazardous. For detailed information please see MSDS.

** The strips of the PT tube can also be evaluated with the device.

Technical specifications QUANTOFIX® Relax REF 91346

Instrument memory

- 200 test results including sample ID

Interface

- User: Touch screen, alphanumeric input, password protection
- Computer: USB and RS 232 interface for connection to PC, PS/2 interface for connection of keyboard or barcode scanner

Power requirements

- 100–240 V AC, automatic
- Battery powered operation (optional) with 6 AA batteries

Dimensions / Weight

- Depth: 20 cm (7.9 inches)
- Width: 16 cm (6.3 inches)
- Height: 7,5 cm (3.0 inches)
- Weight: 710 g (1.90 lb)
(w/o batteries and power supply)

Operating conditions

- Temperature range: 10 °C – 40 °C (50 °F – 104 °F)
- Humidity range: 80 % maximum (non condensing)
- Calibration: automatic, self calibrating

Capacity

- 50 strips per hour



© Ignatius (S. 1) - stock.adobe.com
KATDE100127 QUANTOFIX® Relax en8a

www.mn-net.com

MACHEREY-NAGEL



Management
System
EN ISO 13485:2016
ISO 9001:2015

www.tuv.com
ID 0000056401



MACHEREY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Germany

DE +49 24 21 969-0 info@mn-net.com

CH +41 62 388 55 00 sales-ch@mn-net.com

FR +33 388 68 22 68 sales-fr@mn-net.com

US +1 888 321 62 24 sales-us@mn-net.com

