Phosphatesmo MI

de

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en

Description:

Test strip for the determination of alkaline phosphatase in cow's milk. Phosphatesmo MI is a valuable auxiliary for a quick check of correct pasteurization of milk.

Pack content:

1 aluminum container with 50 test strips
50 PE bags

Limit of sensitivity:

0.5 % raw milk in pasteurized milk

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

Hazard warnings:

This test does not contain hazardous substances that must be labelled.

Instructions for use:

1. Dip test strip briefly into the milk.
2. Shake of excess liquid.
3. In order to prevent the test pad of the test strip from drying out, place the test strip in the enclosed PE bag. Use a new bag for each test to avoid cross-over contamination.
4. Incubate the test strip at 36 °C for 1 h.

Interpretation:

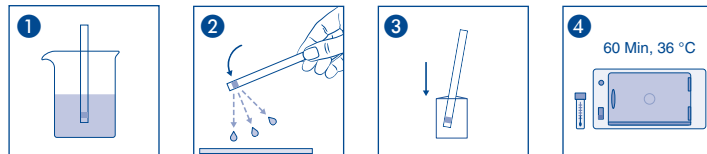
In the presence of alkaline phosphatase the test field turns yellow. The incubation can now be stopped. For control measurement use raw milk and heated milk. A yellow coloration indicates, that raw milk is present or the milk was not sufficiently heated. No coloration indicates that the pasteurization was correctly completed.

Disposal:

Information regarding disposal can be found in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 2–8 °C). If correctly stored, the test strips may be used until the use-by-date printed on the packaging.

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