visocolor® Farbstandard Chlor

Faktor:	1.000
Messwellenlänge (HW = 5 – 12 nm):	530 nm / 540 nm
Messtemperatur:	20 – 25 °C

Methode:

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Tel.: 062 388 55 00 · sales-ch@mn-net.com

visocolor® Color standards Chlorine

Factor: 1.000
 Wavelength (HW = 5 – 12 nm): 530 nm / 540 nm
 Measurement temperature: 20 – 25 °C

Method:

Check solutions to control the photometric accuracy (secondary standards according ISO 9001) for the tests *VISOCOLOR*® Powder Pillows Chlorine free (REF 936220), *VISOCOLOR*® Powder Pillows Chlorine total (REF 936221), *VISOCOLOR*® ECO Chlorine 2, free + total (931015), *VISOCOLOR*® ECO Chlorine 2, free (REF 931016), *VISOCOLOR*® ECO Chlorine 6, free + total (REF 931217) and *VISOCOLOR*® ECO Chlorine 6, free (REF 931219). Suitable for the photometers *NANOCOLOR*® Advance, PF-12 / PF-12^{Plus} and PF-3-series

Content:

- 1 round tube with check solution „Standard 1“
- 1 round tube with check solution „Standard 2“
- 1 round tube with check solution „Standard 3“
- 1 round tube with zero solution "Blank Value"

Procedure:

1. Open the corresponding chlorine-method (e.g. 5 – 17 on the PF-12^{Plus})
2. Perform blank measurement with tube "Blank Value"
3. Insert different test solutions one after another and measure
4. Compare result with certificate

Note:

Tubes must be clean. If necessary, clean them with a dry and clean cloth. Do not use scratched tubes.

Interpretation:

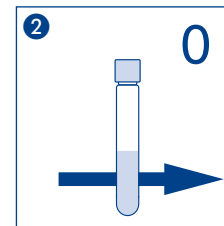
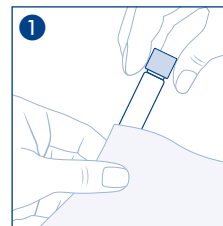
If the measuring values are within the confidence interval, the wavelength settings and measuring geometry of the photometer are in order. If the measuring values are out of the confidence interval, please proceed as follows:

1. Check the power supply
2. Check expiry date of the reagent set; if necessary, use a new pack
3. Check cleanliness of the tubes
4. Check if tubes are scratched
5. Calibrate the photometer and repeat the procedure
6. If necessary, contact your distributor or MACHEREY-NAGEL



Storage:

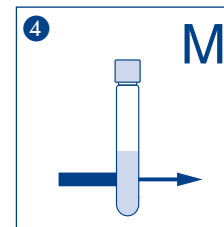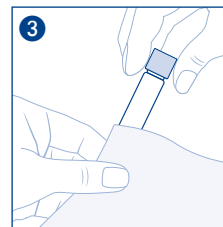Store the check solutions in a cool (15 – 25 °C) place.

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MACHEREY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Alemania