


visocolor® ECO Sauerstoff


Oxygen / Oxygène / Oxígeno

Method(e) / Método	PF-3 Fish	PF-12	PF-12 ^{Plus}	Advance
5881 1-8 mg/L O ₂	530 nm	540 nm	540 nm	540 nm

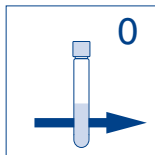


Null mit Probe
Blank with sample
Blanc avec échantillon
En blanco con muestra


Geeignet für Meerwasser
Suitable for seawater
Convient à l'eau de mer
Adecuado para el agua de mar

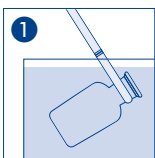

Probe
Sample
Echantillon
Muestra

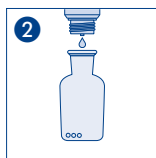

5 ∆ O₂-1

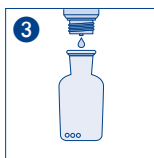

5 ∆ O₂-2



Schütteln
Shake
Agiter
Agitar




Warten
Wait
Attendre
Espere



12 ∆ O₂-3




Schütteln
Shake
Agiter
Agitar

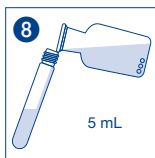

in Rundküvette
into test tube
dans tube
en tubo



Säubern
Clean
Nettoyer
Limpiar

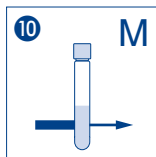

Sofort messen
Measure immediately
Mesure immédiatement
Medir inmediatamente