


visocolor® ECO Chlor 1 (frei)


Chlorine (free) / Chlore (libre) / Cloro (libre)

Method(e) / Método		PF-3 Pool/ Drinking Water	PF-12	PF-12 ^{Plus}	Advance
		530 nm	540 nm	540 nm	530 nm
5351	0.05–2.00 mg/L Cl ₂	■	■	■	■

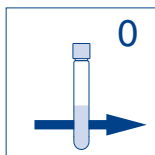


Null mit Probe
Blank with sample
Blanc avec échantillon
En blanco con muestra


Geeignet für Meerwasser
Suitable for seawater
Convient à l'eau de mer
Adecuado para el agua de mar

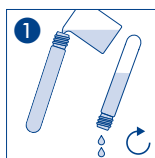


Küvetten spülen
Rinse test tube
Rincer les éprouvettes
Lavar los tubos

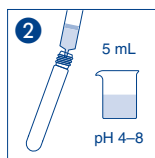


Probe
Sample
Echantillon
Muestra

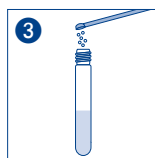

1 Cl₂-1

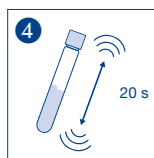


Schütteln
Shake
Agiter
Agitar

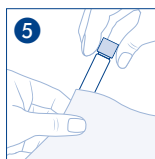

Säubern
Clean
Nettoyer
Limpiar



Warten
Wait
Attendre
Espere

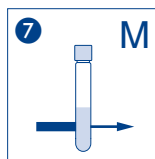


Messung
Measurement
Mesure
Medición

visocolor® ECO Chlor 1 (gesamt)


Chlorine (total) / Chlore (totale) / Cloro (total)

Method(e) / Método	PF-3 Pool / Drinking Water	PF-12	PF-12 ^{Plus}	Advance
5352 0.05–2.00 mg/L Cl ₂	530 nm	540 nm	540 nm	530 nm

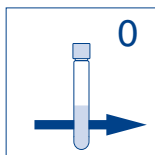


Null mit Probe
Blank with sample
Blanc avec échantillon
En blanco con muestra


Geeignet für Meerwasser
Suitable for seawater
Convient à l'eau de mer
Adecuado para el agua de mar

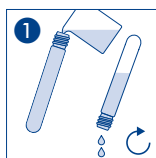


Küvetten spülen
Rinse test tube
Rincer les éprouvettes
Lavar los tubos

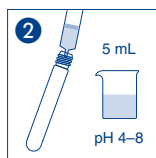


Probe
Sample
Echantillon
Muestra

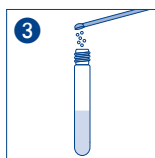


1 Cl_2 -1

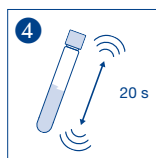


Schütteln
Shake
Agiter
Agitar

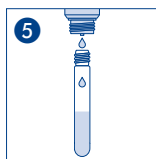

3 Cl_2 -2

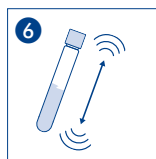


Schütteln
Shake
Agiter
Agitar

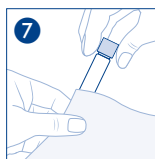

Säubern
Clean
Nettoyer
Limpiar



Warten
Wait
Attendre
Espere

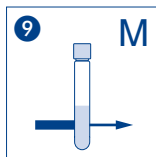


Messung
Measurement
Mesure
Medición

visocolor® ECO Chlor 1 (gebunden)


Chlorine (combined) / Chlore (combiné) / Cloro (ligado)

Method(e) / Método		PF-3 Pool/ Drinking Water 530 nm	PF-12 ^{Plus} 540 nm	Advance 530 nm
5353	0.05–2.00 mg/L Cl ₂	■	■	■


Geeignet für Meerwasser
Suitable for seawater
Convient à l'eau de mer
Adecuado para el agua de mar

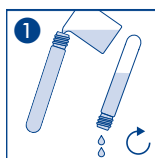


1
Küvetten spülen
Rinse test tube
Rincer les éprouvettes
Lavar los tubos

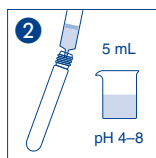


2
Probe
Sample
Echantillon
Muestra

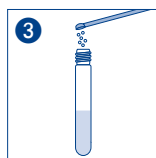

3
1 $\text{Cl}_2\text{-1}$

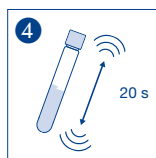


4
Schütteln
Shake
Agiter
Agitar

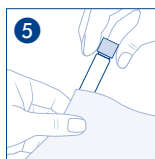


5
Säubern
Clean
Nettoyer
Limpiar




6
Warten
Wait
Attendre
Espere

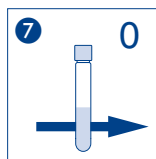


7
Null
Blank
Blanc
Blanco

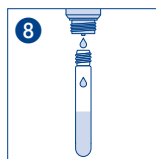

8
3 $\text{Cl}_2\text{-2}$

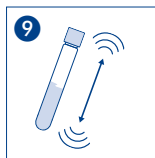


9
Schütteln
Shake
Agiter
Agitar

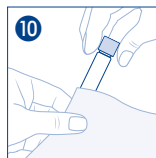

10
Säubern
Clean
Nettoyer
Limpiar



11
Warten
Wait
Attendre
Espere

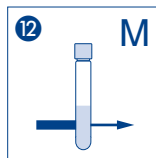

12
Messung
Measurement
Mesure
Medición