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Instructions for using Duotest indicator papers

en

Turn upper half of dispenser anti-clockwise, pull out required strip length, turn clockwise and tear off.

Dip the strip into liquid to be tested.

Compare with color chart on rear of dispenser and read off pH value. Intermediate values may be estimated.

The white center band has been specially impregnated so as to be hydrophobic, thus erroneous readings due to the fusing of indicator dyes, particularly in the alkaline range, are avoided.

Suspensions and viscous solutions should be applied as drops to the indicator paper. For evaluation turn the strip over.

In order to avoid possible errors when weak or nonbuffered liquids are involved, it is advisable to place the strip against a test tube wall. The test tube should then be completely filled. Allow the strip to remain in the liquid for about a minute and compare the color scale with the paper whilst still immersed.

We are continually updating our range of MACHEREY-NAGEL indicator and test papers. Please ask us for full details and do not hesitate to submit to us any special requirements you may have.

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France : MACHEREY-NAGEL SAS · 1, rue Gutenberg – BP135 · 67720 Hoerdt · France
MACHEREY-NAGEL SAS (Société par Actions Simplifiée) au capital de 186600 €