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Rapid Test for determination of protein residues

Contents:

- 1 aluminium container with 60 test strips
- 1 bottle INDIPRO-1
- 1 bottle INDIPRO-2

INDIPRO is a rapid test for determination of protein residues on working surfaces and tools having contact with foodstuffs.

Safety precautions:

Information regarding safety can be found on the box' label and in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

Instructions for use:

1. Add 1 drop of reagent INDIPRO-1 on the test field of the test strip.
 2. Draw test strip with the moistened test field approximately 10 cm under low pressure over the surface to be tested.
 3. Add 1 drop of reagent INDIPRO-2 on the test field.
- A colour change from yellow to green indicates the presence of protein residues. The surface has to be cleaned.

Limit of sensitivity:

50 µg BSA (bovine serum albumin) gives a clear green colouration.

Interference:

Residues of desinfectants can cause also a green colouration of the test field.

Disposal:

Information regarding disposal can be found in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4–30 °C).

If correctly stored, the test kit may be used until the use-by-date printed on the packaging.

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